


## E-Const-Temp Operating Platform Quick Guide


Version 1.0

### Instructions:


1. Take out the instrument and place it on a horizontal and flat surface.
2. Make sure that the power switch of the E-Const-Temp Operating Platform is on the "O" side.
3. Place appropriate metal blocks in the thawing area and low-temperature setting-up area.
4. Connect the power supply and press the power switch to the "I" side, the screen of the instrument will light up and work, and the low-temperature setting-up area will start to cool down. At the same time, the thawing area began to heat up.
5. After the low-temperature setting-up area reaches the working temperature, the screen will display "Ready", and the experiment can be carried out.
6. At the end of the experiment, press the power switch to the "O" side and pull out the power plug from the socket.

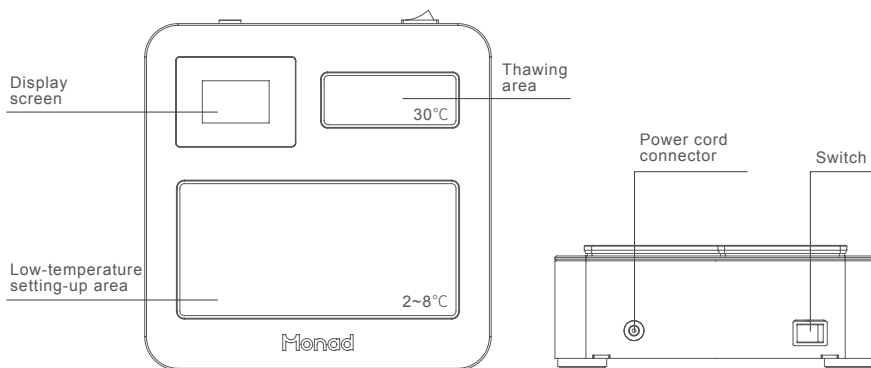

### 1. Screen display instructions

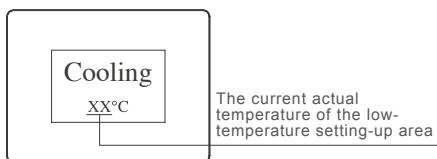

Preparation phase of low-temperature setting-up area

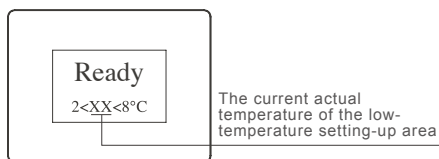

Working status of low-temperature setting-up area